


Mouse and Rat Restraints

Traditional Triangle Shape

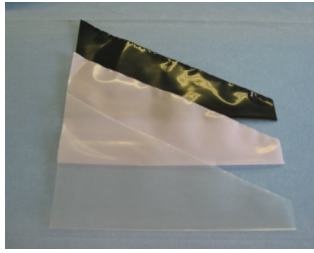


*Includes instructions
& twist ties*

Size	Animal	Clear	Black	Rose
X-Small <i>mic</i>	< 20 g	1800C	xx	xx
Small	≤ 120 g	1801C	1801B	1801R
Medium	≤ 225 g	1802C	1802B	1802R
Large	≤ 450 g	1803C	1803B	1803R

● 50 restraints/pack, custom sizes available ●

Humane
Safe
Disposable

Body-Contour™ Cone Shape

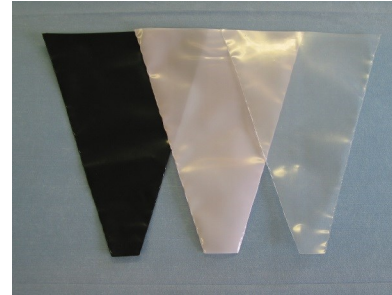


Size	Animal	Clear	Black	Rose
Small	≤ 120 g	1801CV	1801BV	1801RV
Medium	≤ 225 g	1802CV	1802BV	1802RV
Large	≤ 450 g	1803CV	1803BV	1803RV


● 40 restraints/pack, custom sizes available ●

In Low Tail Injection Platform

Cat #: IL300

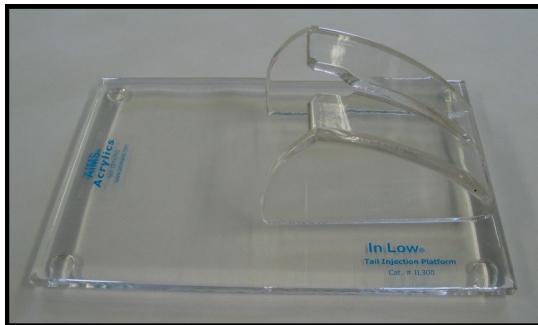


- ◆ Quick Restraint
- ◆ Low Stress
- ◆ Easy to Clean

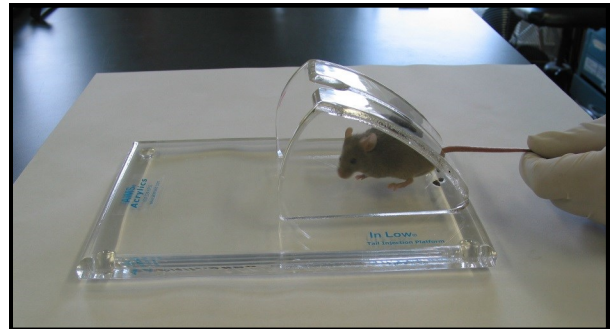


In Low Tail Cutting Platform

Cat #: IL200

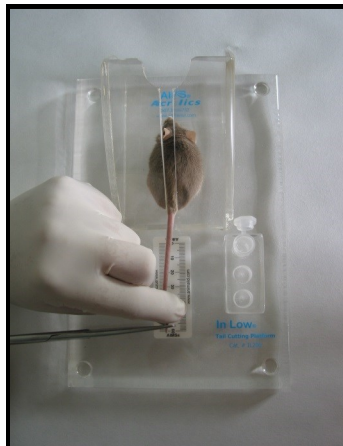

- ◆ Quick Restraint
- ◆ Tube Holder
- ◆ Ruler for Accurate Snips
- ◆ Easy to Clean

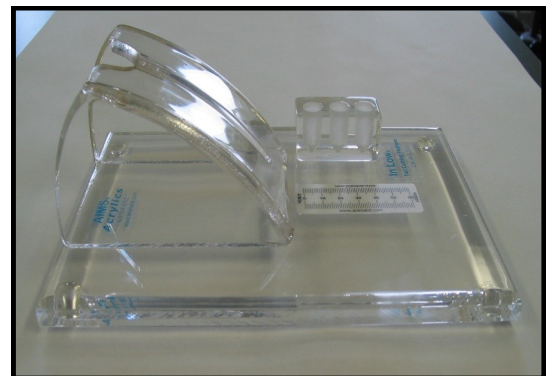


Order using Visa, MasterCard, American Express or an Official PO

AIMS Rodent Restraint Bags



Step 1

Hold large end of bag open and allow rat to enter.

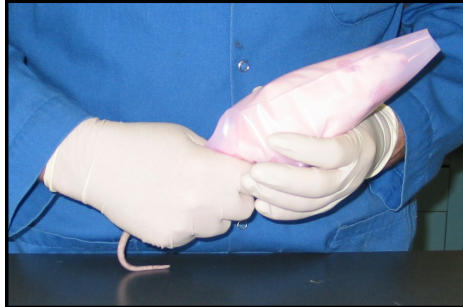

Step 2

As rat enters, keep narrow end of **bag & animal pointed up**. *Animals prefer to run upwards.*



Step 3

Carefully twist plastic bag around base of tail and use twist tie to secure plastic to tail.

Technical Notes:

- Do not leave animal in bag for extended periods of time. Animal body heat will increase when restrained in a plastic bag.
- Placing a mark at the base of the tail (on top) before placing animal in bag will help with tail orientation when animal is fully immobilized.
- Black/Rose colored restraint bags reduce light and provide a more calming environment for animals.
- Body-Contour cone shape bags provide easier animal capture and increased animal comfort.



Voice 607.324.6752 • www.animalid.com
aims@animalid.com • Fax 607.324.6753

All text and graphic are © 2021 by Animal Identification
and Marking Systems, Inc. All rights reserved.